


食品分析

3-1

37 种顺式/反式脂肪酸甲酯的分析
(以 50m 长度的柱子解决其它公司 100m 长度柱子的分离)

组份: 37 种顺式/反式脂肪酸甲酯

色谱柱: MEGA-10 50m x 0.25mm 0.25um df
订货号: C-10-025-025-50

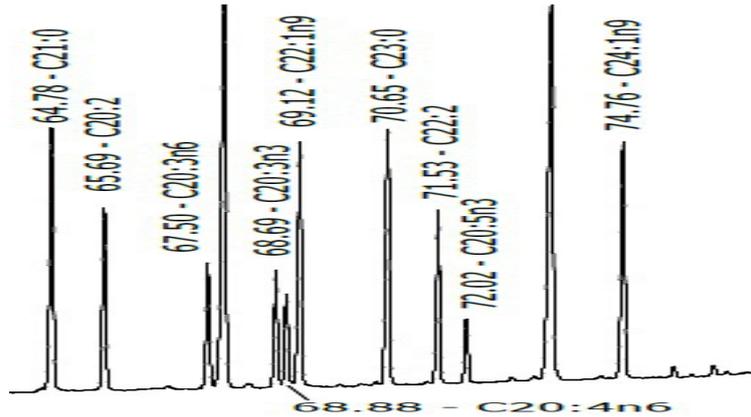

MEGA
CAPILLARY COLUMNS
LABORATORY

Conditions:
Injection: Split 250°C, 1µL, 1:100 Split Ratio.
Sample: 37 FAMES Mix (Supelco cat. #: 47885-U).
Detector: FID 255°C.
Oven Temp.: 60°C (6min), 2°C/min, 230°C.
Carrier Gas: Hydrogen, 100 kPa.

FAMES Mix
(*cis/trans* Isomers) on MEGA-10
Column: **MEGA-10 - 0.25mm, 0.25µm, 50m**
Catalog Code: C-10-025-025-50

Zoom In to see the chromatogram in high definition and better appreciate the separation!

since 1980
MEGA s.n.c. - Capillary Columns Laboratory - Via Plinio, 29 - 20025 Legnano (MI) - Italy - Tel./fax: (+39) 0331.54.79.24 - info@mega.mi.it
www.mega.mi.it

